

DIRECTIONS FOR FITTING BEAM COVER ON #29 BALANCE

1. Remove beam and bows
2. Remove two bearing screws that hold the lift rods in the arms.
3. In the top of the column there are two screws that hold the cap or stone piece on the column. Remove these screws and lift cap off the column. The arms being fastened on the cap, they will come off also.

CAUTION There are four screws, two on each side that hold the arms on the cap. These screws must not be removed under any conditions.

4. Cap and arms having been removed in one unit, the bottom of the beam cover should be slipped over the top of the column and over the lift rods, until the brass cover support rests on the small shoulder on the column provided for that purpose.

Care should be taken to have the brass cover support on the upper side of the board nearest the top of the column.

5. Replace cap and arms and screw firmly in place. Replace lift rods in the arms.
6. Replace beam and bows and place cover on the board.